



SMOBio
Small Bio, Smart Tool

www.smobio.com

Product Information

ExcelBand™

**1 KB Plus (0.1-10 kb) DNA Ladder,
Ready to Use**

DM3200 87 µg / 500 µl

Size: 500 µl

5 µl per well for 100 gel applications

Storage

Room temperature	≥ 6 months
4°C	≥ 12 months
-20°C	≥ 24 months

Description


The DM3200 ExcelBand™ 1 KB Plus (0.1-10 kb) DNA Ladder is a ready-to-use DNA ladder, which is pre-mixed with loading dye for direct gel loading. The DNA Ladder DM3200 is composed of 19 individual DNA fragments: 10k, 8k, 6k, 5k, 4k, **3k**, 2.5k, 2k, **1.5k**, 1k, 900, 800, 700, 600, **500**, 400, 300, 200, and 100 bp derived from a mixture of PCR products and specifically digested plasmid DNA. This product contains three enhanced bands (3 kb, 1.5 kb and 500 bp) for easy reference. In addition, three tracking dyes, Xylene cyanol FF, Bromophenol blue and Orange G which mimic the migration of 4,000 bp, 500 bp and 50 bp dsDNA during electrophoresis are added for real time monitoring.

Buffering

10 mM Tris-HCl (pH 8.0), 10 mM Na-EDTA, 0.025% Orange G, 0.005% Bromophenol blue 0.005%, Xylene cyanol FF and 10% Glycerol.


Quality Control

1. 19 defined bands are observed during electrophoresis in agarose gel.
2. No nuclease activity is detected by a direct nuclease activity assay.

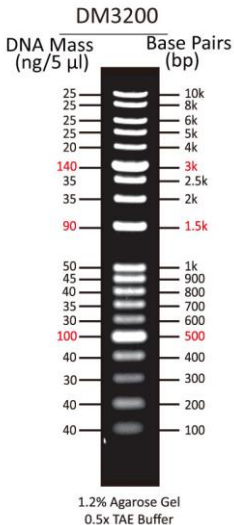

Note: This DNA Ladder was not designed for precise quantification of DNA mass, but can be used for approximating the mass of DNA by intensity comparison.

ExcelBand™ 1 KB Plus (0.1-10 kb) DNA Ladder

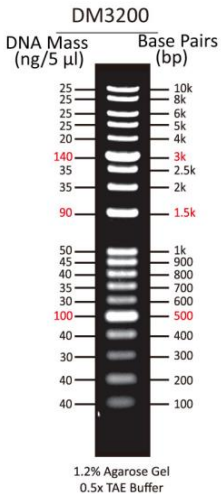

The ladder is supplied in a gel loading buffer and is ready to use.
Do not heat or dilute before loading.

Other information

SMOBIO Technology, Inc. claims all warranties with respect to this document, expressed or implied, including but not limited to those of merchantability or fitness for a particular purpose. In no event shall SMOBIO Technology, Inc. be liable, whether in contract, tort, warranty, or under any statute or any other basis for special, incidental, indirect, punitive, multiple or consequential damages in connection with or arising from this document, including but not limited to the use thereof.

Caution: Not intended for human or animal diagnostic or therapeutic uses.

Related Products

DM1100	ExcelBand 50 bp DNA Ladder, 500 μ l
DM2100	ExcelBand 100 bp DNA Ladder, 500 μ l
DM2300	ExcelBand 100 bp+3K DNA Ladder, 500 μ l
DM3100	ExcelBand 1 KB (0.25-10 kb) DNA Ladder, 500 μ l
DM4100	ExcelBand XL 25 kb DNA Ladder, Broad Range (up to 25 kb), 500 μ l
DM5100	ExcelBand Super Range DNA Ladder, (50 bp- 25 kb), 500 μ l
DL1000	ExcelDye DNA Loading Dye, Orange, 5 ml \times 2
DL2000	ExcelDye DNA Loading Dye, Green, 5 ml \times 2
DL3000	ExcelDye DNA Loading Dye, Blue, 5 ml \times 2
DL4000	ExcelDye DNA Loading Dye, Tri-color, 5 ml \times 2
DL5000	FluoroDye DNA Fluorescent Loading Dye (Green, 6X), 1 ml
DL5001	FluoroDye DNA Fluorescent Loading Dye (Green, 6X), 1 ml \times 5
DS1000	FluoroStain DNA Fluorescent Staining Dye (Green, 10,000X), 500 μ l
NS1000	FluoroVue Nucleic Acid Gel Stain (10,000X), 500 μ l
VE0100	B-BOX TM Blue Light LED epi-illuminator



VE0100 B-BOX™ Blue Light LED epi-illuminator

For research use only
2016 ver. 1.3.0