



**SMOBIO**

Small Bio, Smart Tool

[www.smobio.com](http://www.smobio.com)

## Product Information

### SMO-HiFi™ DNA Polymerase

#### TF1000 100 units

SMO-HiFi™ DNA Polymerase (1 U/μl)	100 μl
10X HiFi™ Buffer	600 μl
25 mM MgSO <sub>4</sub>	500 μl
dNTPs Mix (2 mM each)	600 μl
DMSO	600 μl

## Storage

-20°C for 24 months

## Applications

- High fidelity PCR
- Generates blunt end amplicons for cloning with GetClone™ PCR cloning vector

## **Description**

The SMO-HiFi™ DNA Polymerase is a new genetically modified, recombinant DNA polymerase with 70 times higher fidelity than *Taq* DNA polymerase during amplification, as well as very high elongation rate. Being highly thermostable, SMO-HiFi™ DNA Polymerase can remain viable even after being subjected to boiling for 2 minutes. The SMO-HiFi™ DNA Polymerase is also designed to operate in much lower  $Mg^{2+}$  concentration as compared to other DNA polymerase products.

## Features

- 5'→3' DNA polymerase activity
- 3'→5' exonuclease (proofreading) activity
- High reaction rate: 10 seconds/kb
- High fidelity: 70 times higher than *Taq* polymerase
- Generates blunt end amplicon
- Thermo-stable for more than 10 hrs at 95°C.

## Storage Buffer

50 mM Tris-HCl (pH 8.0), 50 mM KCl, 0.1 mM EDTA, 1 mM DTT, stabilizer, 50% (v/v) glycerol

## Unit Definition

One unit is defined as the amount of enzyme that will incorporate 10 nmol of dNTP into acid-insoluble material in 30 minutes at 74°C.

## Recommended PCR Condition

Template	1 – 150 ng
Forward primer	0.1 – 0.5 $\mu$ M
Reverse primer	0.1 – 0.5 $\mu$ M
10X HiFi™ Buffer	5 $\mu$ l
MgSO <sub>4</sub> (25 mM)	2 $\mu$ l
dNTPs (2 mM each)	5 $\mu$ l
SMO-HiFi™ DNA Polymerase	1 $\mu$ l (1 unit)
DMSO	5 $\mu$ l
H <sub>2</sub> O	to 50 $\mu$ l
<hr/>	
Total volume	50 $\mu$ l

## Recommended PCR Program

94°C	2 min	} 25 ~ 40 cycles
94°C	15 sec	
50~68°C*	30 sec	
68°C	30 sec/kb	
68°C	1 min	

\*Optimal PCR condition varies according to primers' thermodynamic properties.

## **Quality Control**

### **Functional Testing**

SMO-HiFi™ DNA Polymerase is tested for performance in the polymerase chain reaction (PCR) using 1 unit of enzyme to amplify a 665 bp target from 1 pg of tested plasmid DNA. The resulting PCR product is visualized as a single band on an ethidium bromide-stained agarose gel.

### **Nuclease Assay**

No contaminating endonuclease or exonuclease activity was detected using pUC19 incubated with SMO-HiFi™ DNA Polymerase for 4 hours at 37°C.

### **Residual Nucleotides Assay**

No contaminating residual nucleotides were detected from purified SMO-HiFi™ DNA Polymerase by PCR assay.

## **Other Information**

SMOBIO Technology, Inc. claims all warranties with respect to this document, expressed or implied, including but not limited to those of merchantability or fitness for a particular purpose. In no event shall SMOBIO Technology, Inc. be liable, whether in contract, tort, warranty, or under any statute or any other basis for special, incidental, indirect, punitive, multiple or consequential damages in connection with or arising from this document, including but not limited to the use thereof.

Caution: Not intended for human or animal diagnostic or therapeutic uses.

## Related Products

CK1000	Champion E. coli Transformation Kit
CV1000	GetClone PCR Cloning Vector, 20 RXN
CV1100	GetClone PCR Cloning Vector II, 20 RXN
DM2100	ExcelBand 100 bp DNA Ladder, 500 $\mu$ l
DM2300	ExcelBand 100 bp+3K DNA Ladder, 500 $\mu$ l
DM3100	ExcelBand 1 KB (0.25-10 kb) DNA Ladder, 500 $\mu$ l
DL5000	FluoroDye DNA Fluorescent Loading Dye (Green, 6 $\times$ ), 1 ml
NS1000	FluoroVue Nucleic Acid Gel Stain (10,000X), 500 $\mu$ l
TP1000	ExcelTaq DNA Polymerase, 500 U
TP1100	ExcelTaq 5 $\times$ PCR Master Mix, 200 RXN
TP1200	ExcelTaq 5 $\times$ PCR Master Dye Mix, 200 RXN
TP1260	ExcelTaq 5 $\times$ Fluorescent PCR Master Mix, 200 RXN
TP2000	ExcelTaq Blood Direct DNA Polymerase, 500 U
TP5000	ExcelTaq Hot Start II DNA Polymerase, 500 U
TF1100	SMO-HiFi 2X PCR Master Mix, 50 RXN
TF2000	Q-HiFi DNA Polymerase, 1 U/ $\mu$ l, 100 U
TF3000	G-HiFi DNA Polymerase, 1 U/ $\mu$ l, 100 U
VE0100	B-BOX Blue Light LED epi-illuminator, AC 100-240V, 50/60Hz



B-BOX™ Blue Light LED epi-illuminator

For research use only  
2017 ver. 2.1.0